

## Gamma Wipe® - VP

Sterile Sealed Edge Cleanroom Laundered 100% Polyester Knit wiper

Gamma Wipe® VP is an ultrasonically sealed edge cleanroom laundered wiper recommended for ISO Class 3 and above environments composed of 100% continuous filament polyester knit fabric. Each package is gamma irradiated sterile to a 10<sup>-6</sup> Sterility Assurance Level. This combination of properties provides Gamma Wipe® VP with the highest level of cleanliness, abrasion resistance and chemical compatibility required for cleaning critical sterile processing environments.



### Key Attributes

- 100% continuous filament polyester knit
- Ultrasonically sealed edge for reduced fiber contamination
- Laundered and packaged in Berkshire's ISO Class 4 cleanroom with Lot Number on each pack
- Gamma irradiated and sterile validated to a 10<sup>-6</sup> Sterility Assurance Level per AAMI guidelines
- Linear tear packaging
- Certificate of Sterility with each case and available on the web by lot number 24/7. Details: Date of Manufacture, Radiation Dosage, Lot Information, Meets cGMP requirements for traceability

### Benefits

- Critically low particles, fibers, ions and extractables
- High abrasion resistance
- High Absorbency
- Chemically compatible with common cleaning and disinfecting solutions including isopropyl alcohol, phenols, bleaches and quarternary ammonium compounds
- Light weight, high absorbency material design reduces actual cost in use compared to heavier weight products

### Environmental

- Light weight, high absorbency design reduces landfill waste impact when compared to more traditional heavier weight designs
- Complies with RoHS

### Applications

- Designed for use in ISO Class 3 and higher sterile environments where microbial exclusion is demanded
- Designed for cleaning sterile surfaces and equipment during aseptic processing and USP <797> applications
- Excellent for cleaning sterile laminar air flow units, barrier isolators and bio-containment units

### Pre-Wetted Option

The same wiper material can be provided in pre-wetted formats for reduced VOC emissions, increased convenience, increased productivity, improved solvent control and cleaning protocol repeatability and reduced costs. Pre-wetted options include WFI.

### Validated Sterile

ANSI/AAMI/ISO procedures have been followed for verification of irradiation dose with quarterly audits being performed to monitor bioburden levels and dose. The Sterility Assurance Level (SAL) or the probability of a viable organism being present on a product unit after sterilization is a SAL of 10<sup>-6</sup> (which is the chance of 1 non-sterile unit in 1,000,000).

### Other Sterile wipers

- Gamma Wipe® 300
- Gamma Wipe SuperSorb®
- Gamma Wipe®120
- Gamma Wipe® 67
- Gamma Wipe® CapSure™-VP

[www.berkshire.com](http://www.berkshire.com)

Contact: Tel 1 800 242 7000 / 1 413 528 2602  
 info@berkshire.com

America	Tel 1 413 528 2602	Fax 1 413 528 2614	info@berkshire.com
Europe	Tel 44 (0) 870 757 2877	Fax 44 (0) 870 757 2878	enquiries@berkshire.uk.com
SE Asia	Tel 65 6252 4313	Fax 65 6252 4312	enquiries@berkshire.com.sg
Japan	Tel 81 3 5827 2380	Fax 81 3 5827 2382	master@berkshire.co.jp

## Technical Data:

Attribute		Units	Value	Test Method
<b>Basis Weight</b>		g/m <sup>2</sup>	<b>130</b>	TAPPI T-410
<b>Caliper</b>		um	<b>442</b>	TAPPI T-411
<b>Fibers</b>	>100um	fibers/cm <sup>2</sup>	<b>0.12</b>	IEST-RP-CC004.3, Sec 6.1.3 / Sec 6.2.2
<b>Particles</b>	>0.5um	x10 <sup>3</sup> /cm <sup>2</sup>	<b>0.31</b>	IEST-RP-CC004.3, Sec 6.1.3 / Sec 6.2.1
<b>Sorbency</b>	Capacity	mL/m <sup>2</sup>	<b>448</b>	IEST-RP-CC004.3, Sec 8.1 modified / Sec 8.2 modi-
	Efficiency	mL/g	<b>3.4</b>	
	Rate	seconds	<b>1</b>	
<b>Non-Volatile Residue</b>	DI Water	g/m <sup>2</sup>	<b>0.0019</b>	IEST-RP-CC004.3, Sec 7.1.2
	IPA	g/m <sup>2</sup>	<b>0.0054</b>	
<b>Ions</b>	Na <sup>+</sup>	ppm	<b>0.17</b>	IEST-RP-CC004.3, Sec 7.2.2
	K <sup>+</sup>	ppm	<b>0.024</b>	
	Ca <sup>++</sup>	ppm	<b>0.053</b>	
	Mg <sup>++</sup>	ppm	<b>0.011</b>	
	Cl <sup>-</sup>	ppm	<b>0.032</b>	

### Notes:

- All values are statistical averages
- Due to differences in test methods applied and equipment utilized by different wiper manufacturers, valid product comparisons may only be obtained through side-by-side testing in the same test facility, under similar conditions
- Third party testing can be performed upon request

## Order Information:

Product	Number	Size	Shts/pk	Pks/Inner bag	Inner bags/cs	Pks/cs	Style
Gamma Wipe <sup>®</sup> VP	GWVPST0940	9x9"(23x23cm)	20	5	8	40	Header Bag
Gamma Wipe <sup>®</sup> VP	GWVPST25	12x12"(30x30cm)	20	5	5	25	Header Bag

## Other Berkshire Products




Wipers



Glove Liners



Mop Systems

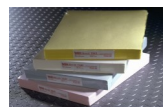

Documentation Systems



Face Masks



Swabs