Analyzer for Water Quality & Food

BOD Meter
COD Meter
MLSS Meter
AMMONIA Meter
H₂O₂ Meter
FRESHNESS METER

BOD analyzer with Biosensor

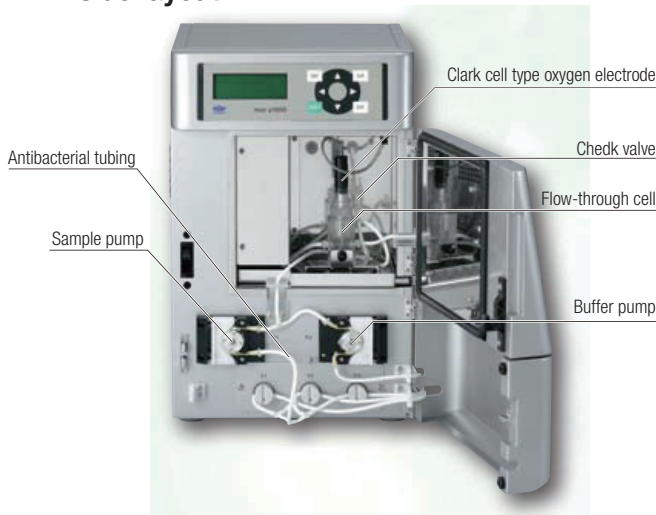
Quick BOD α 1000

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Inside layout

Easily viewable display(20 digits x 4 lines, back-light LCD). Simple operating keys.

Specifications

측정방법	Biosensor
DO 전극	Clark cell type oxygen electrode
측정범위	2~50mg/L (basic measurement range based on trichosporon cutaneum membrane)
측정시간	Approx. 60 minutes per measurement
측정 가능한 샘플	Liquid sample containing soluble organic
캘리브레이션	1 point calibration by glucose glutamic solution
재현성	±5% (in case of 20~50ma/L F.S.setting, using trichosporon cutaneum membrane and stable environment temperature)
온도조절	20 to 40 (adjustable) Automatic warming/cooling PI control function with peltiert device
외부출력	Print-out output BOD value hold output (DC 0~5V) Probe output (DC 0~5V)
공급전원	100VAC, 50/60Hz, 3A
크기	206(W) x 320(D) x 409(H) mm
중량	Approx. 16kg, not including reagents
약세서리(옵션)	Microbe membrane (trichosporon cutaneum) Pipettes or measuring apparatus

Instrument

Product	SET configurations	Quantity
BOD 1000 Set	<ul style="list-style-type: none"> • Printer • Microbial immobilization kit • BOD seeds • DO Probe Kit • 0.5M Phosphate buffer solution • 5000 mg/L BOD standard solution 	1 set

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Instrument


Order No.	Descriptions	Model	Remarks
C00018301	COM Meter	HC-607	Acid method
C00018302-2	COM Meter	HC-607 with printer	Acid method solution A and B are not included

Consumables

Order No.	Descriptions	Order No.	Descriptions
C00060601	COD Reagent A solution	C00042203	PVC rod
C00060602	COD Reagent B solution	C00042119	Heat-control plate
C00060611	COD Reagent set • 1 Reagent A solution • 2 Reagent B solution	C00042118	Special heater
C00102205	Watch glass Ø 60	C00042120	Special heater with heat-control plate
C00042101	Special tall beaker, 100mL	C00023106	Electrolysis electrode
C00042103	Safety bands	C00023107	Indicator electrode
C00042202	Emery (100g)	C00041411	Printer paper

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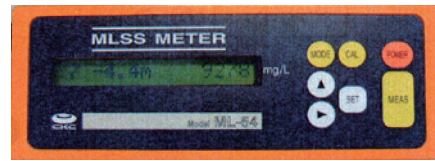Instrument

Order No.	Descriptions	Model	Remarks
C00005410	Portable MLSS meter set	ML-54	Measurement range : 0~15,000 mg/L
C00005411	Portable MLSS meter set with printer	ML-54	Measurement range : 0~15,000 mg/L
C00005412	Portable MLSS meter set	ML-54	Measurement range : 0~15,000 mg/L
C00005414	Portable MLSS meter set with printer	ML-54	Measurement range : 0~15,000 mg/L
C00031120	MLSS sensor, ML-54		Sensor cable 10m
C00031121	MLSS sensor, ML-54		Sensor cable 15m
C00042225	Printer for MS-54	SD1-31SJ	Interface cable (straight)
C00051112	Carrying case, ML-54		Aluminium

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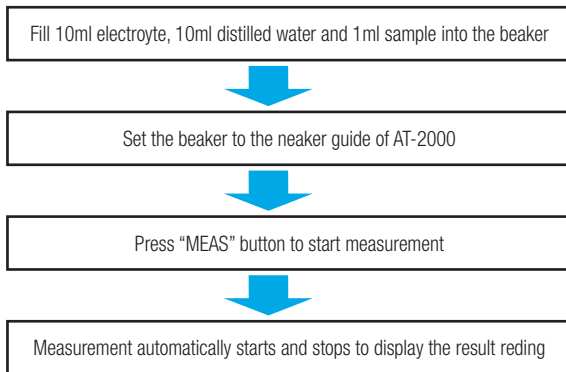
Easy steps for the sample measurement (for the case of the range 0-20mg/L)

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
Result data of the standard solution measurement

Concentrarion(mg/L)	Instrument reading(mg/L)
0.2	0.21
1	1.02
2	2.02
10	9.94
20	19.9

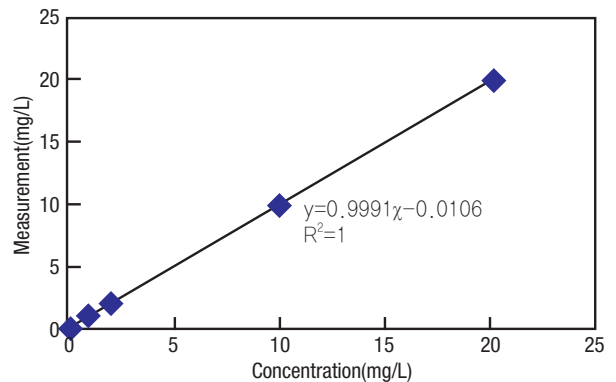Average of 3 repetitions is shown for each concentration.
Average CV was 0.78%

Instrument

Order No.	Descriptions	Model	Remarks
C00017211	Quick Ammonia	AT-2000	" • Analyzer • Eletrolyte (2) • Printer

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Instrument

Order No.	Descriptions	Model	Remarks
C00250001	H2O2 Meter, Super Oritector	M-5	• Meter • Reagents
C00250401	Printer for M-5		• Printer connection cable included
C00250201	Constant-temperature Bath for M-5		
C00051112	Carrying case, ML-54		

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Instrument

Order No.	Descriptions	Model	Remarks
C00250301	Freshness Meter	KV-202	• Meter • Reagents • Printer • Constant - temperature Bath
C00250401	Printer for KV-202		• Printer connection cable included
C00250201	Constant-temperature Bath for KV-202		

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