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\*Revodix Inc  
#308 CRESCO Bldg, 114, Kwangjang-dong,  
Kwangjin-ku, Seoul, Korea. 143-802

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Note: The step resolution sated is the minimum step resolution achieved by a pump.

### **Fusion Touch 100**

Syringe size	0.5 microliter to 60 milliliter
Voltage operating range	115 V~240V, 0.25 A, 55-60 Hz
Drive mechanism	Stepper motor
Step resolution (advance per microstep)	0.198 microns
Flow rate range	0.001 $\mu$ l/hr (10 $\mu$ l) to 90 ml/min
Nominal linear force	35 lbs
Dimensions	9.5x 6.5 x 4 inch
Weight	7.0 lbs
Temperature range	10°C ~ 50°C
Humidity	20% - 80% RH
RS232 Connector	D9 Sub-connector
TTL Connector	USB-B Receptacle

### **Fusion 200 Touch**

Syringe size	0.5 microliter to 60 milliliter
Voltage operating range	115 V~240V, 0.25 A, 55-60 Hz
Drive mechanism	Stepper motor
Step resolution (advance per microstep)	0.046 microns
Flow rate range	0.001 $\mu$ l/hr (10 $\mu$ l) to 45 ml/min
Nominal linear force	50 lbs
Dimensions	9.5x 6.5 x 4.5 inch
Weight	7.0 lbs
Temperature range	10°C ~ 50°C
Humidity	20% - 80% RH
RS232 Connector	D9 Sub-connector
TTL Connector	USB-B Receptacle

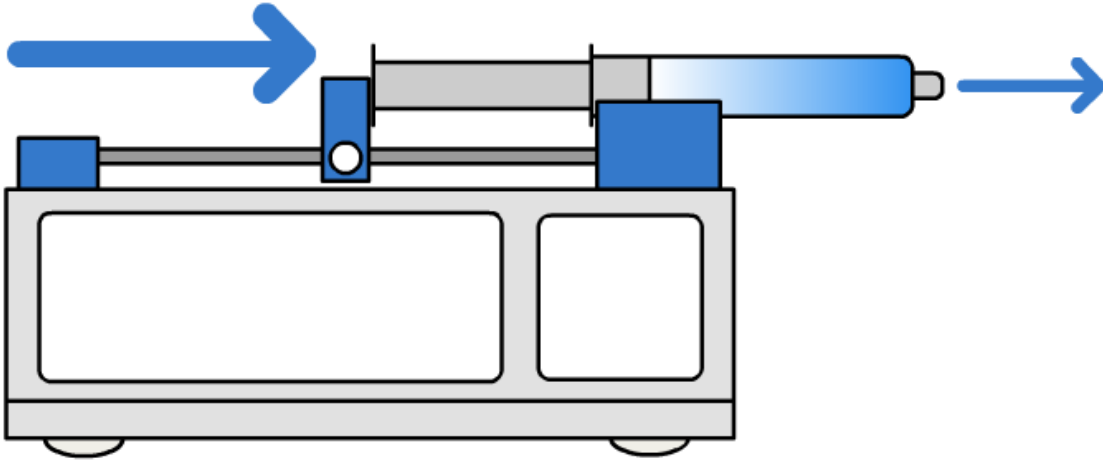
### **Fusion 400 Touch**

Syringe size	0.5 microliter to 10 milliliter
Voltage operating range	115 V~240V, 0.25 A, 55-60 Hz
Drive mechanism	Stepper motor
Step resolution (advance per microstep)	0.016 microns
Flow rate range	10pl/min (0.5 $\mu$ l) to 10 ml/min
Nominal linear force	50 lbs
Dimensions	9.5x 6.5 x 5.5 inch
Weight	7.5 lbs
Temperature range	10°C ~ 50°C
Humidity	20% - 80% RH
RS232 Connector	D9 Sub-connector
TTL Connector	USB-B Receptacle

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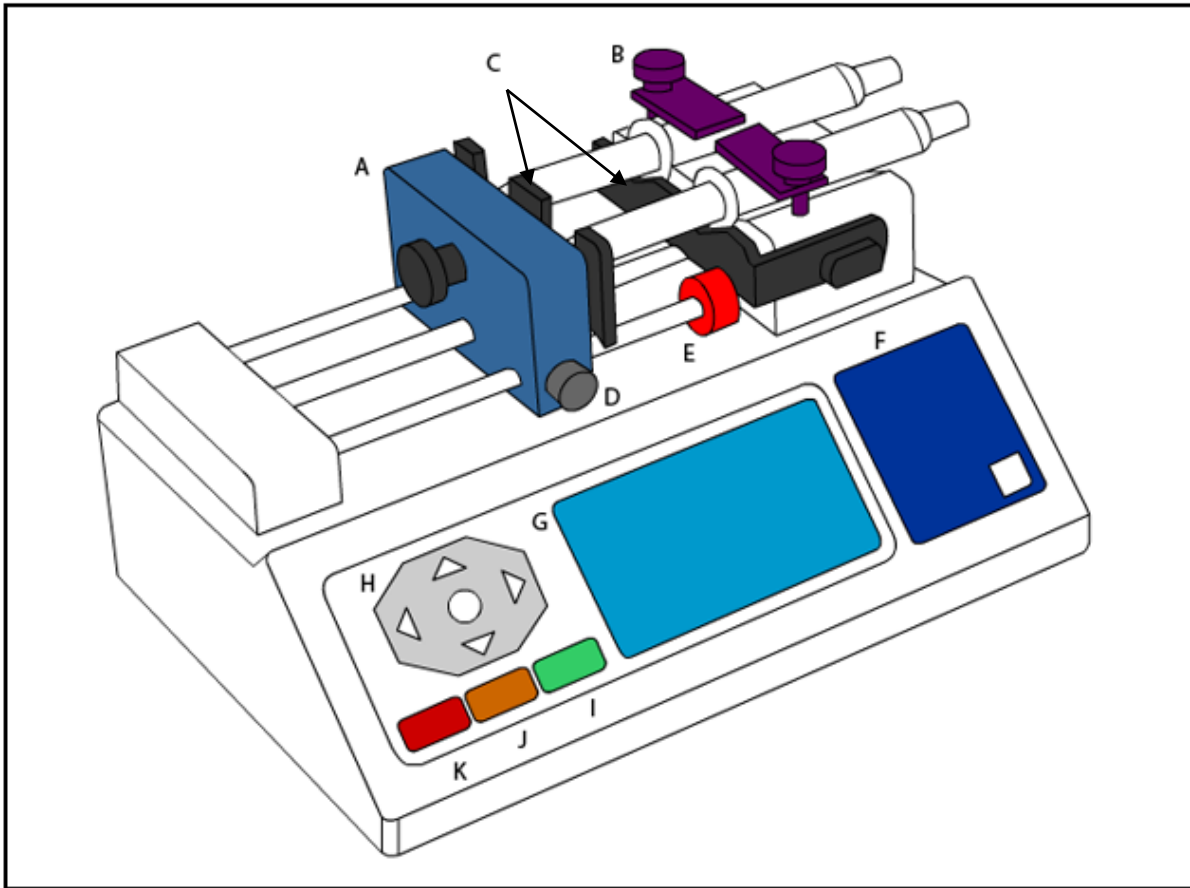
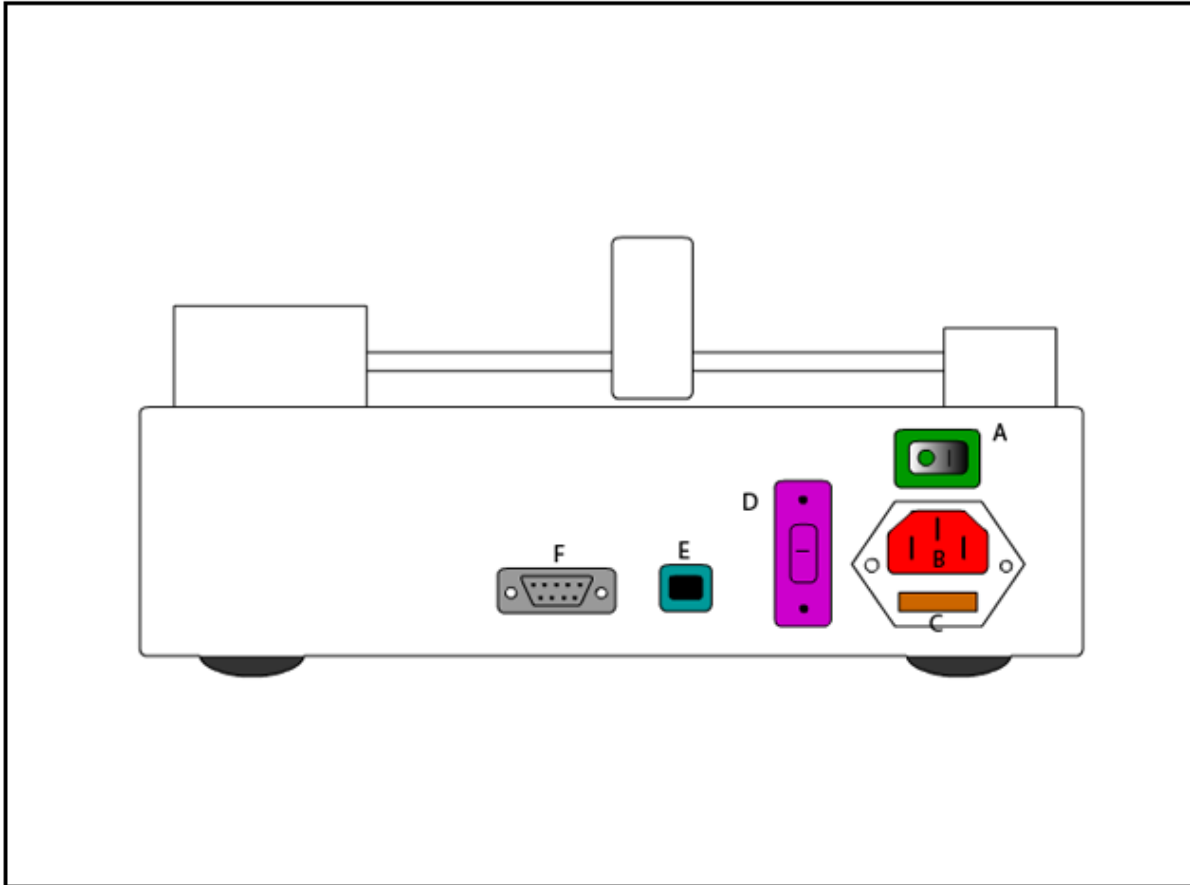
#### **Multi-Step (Optional Multi-step Firmware Upgrade):**

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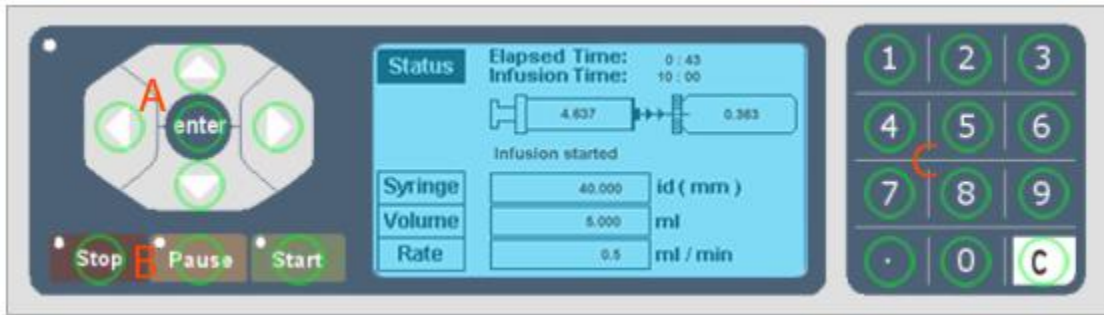
#### **Looping (Optional Multi-step Firmware Upgrade):**

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#### **Ramping Rate (Optional Multi-step Firmware Upgrade):**

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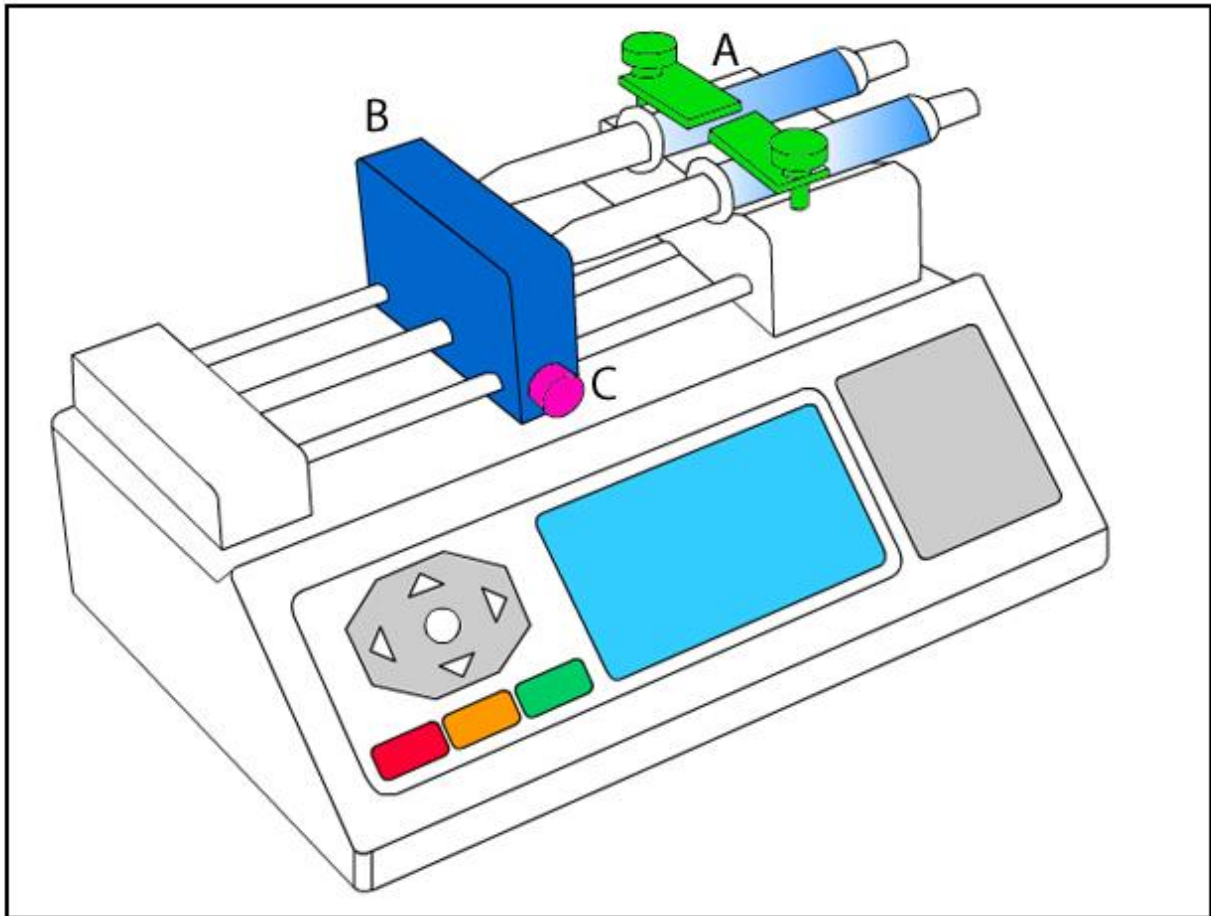
5

Chemyx Inc. <http://www.chemyx.com>

All copyrightable text and graphics, arrangement, and presentation of all materials and overall design of this pump interface are ©2010 Chemyx Inc. All rights reserved

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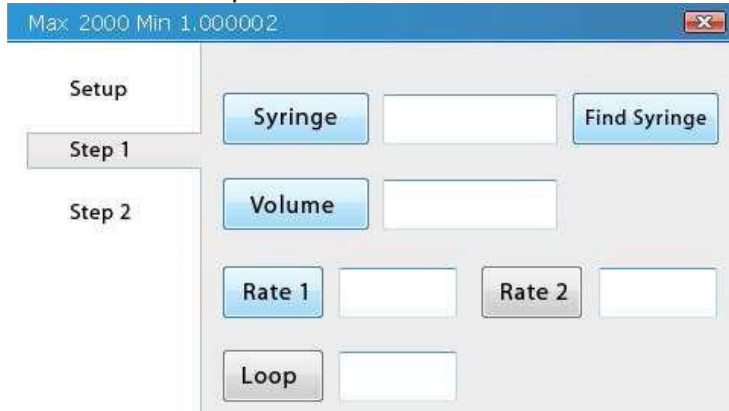
## Flow Profile Setup

The image shows a software window titled "Flow Profile Setup" with a blue title bar containing "Max 2000 Min 1.000002" and a close button. On the left, there is a vertical menu with "Setup", "Step 1", and "Step 2". The main area contains several input fields and buttons: a "Syringe" button next to an empty text box, a "Find Syringe" button, a "Volume" button next to an empty text box, "Rate 1" and "Rate 2" buttons each next to an empty text box, and a "Loop" button next to an empty text box.

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### **Approved Lubricants:**

Engine oil or Motor oil – any type  
Silicone oil  
Machine oil  
Tooling oil or “WD-40”

Contact Chemyx if you wish to use other lubricants.

### **Signs that your system is not adequately lubricated:**

- 1) Grinding sounds coming from the lead screw.
- 2) Locking nut is decoupling under load before stall
- 3) Slow decrease of max pushing force.

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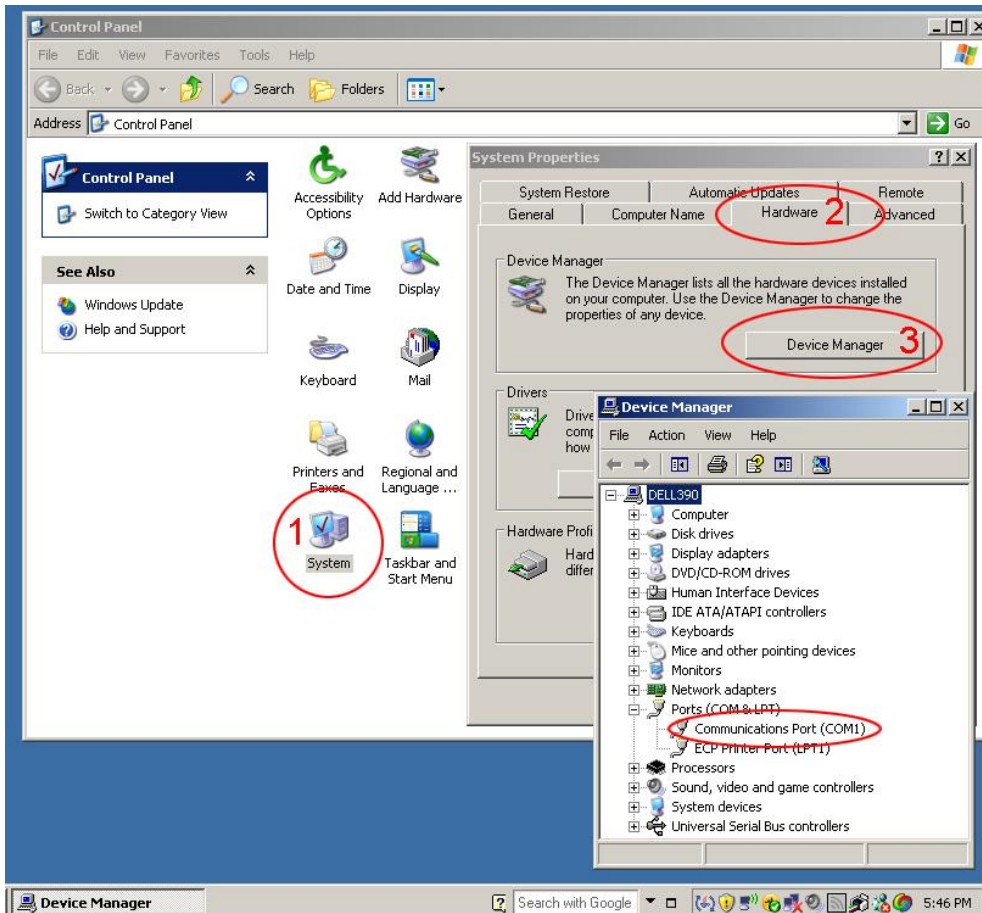
DB9 Serial Cable Male to Female – Strait though configuration.

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Baud Rate - 9600, Data Bits - 8, Parity - none, Stop bits - 1, Flow control – none.

모든 프로그래밍 패키지 - LabView, LabWindows 와 Visual Studio will allow you to program comport settings dynamically in program.

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<u>Protocols:</u>	<u>Resulting Action by Pump</u>
help	- Show help information
stop	- Pump Stop
pause	- Pump Pause
start	- Pump Run
status	- Pump Return Status
set units [xxx]	- Pump Set units
set diameter [x.x]	- Pump Set diameter
set rate [x.x]	- Pump Set rate
set time [xxx]	- Pump Set time@volume
set volume [x.x]	- Pump Set volume
set delay [xxx]	- Pump Set delay@start
save setting	- Pump parameter saved
read limit parameter	- Pump Returns Limit Parameter
view parameter	- Pump parameter settings
restart	- Pump restart

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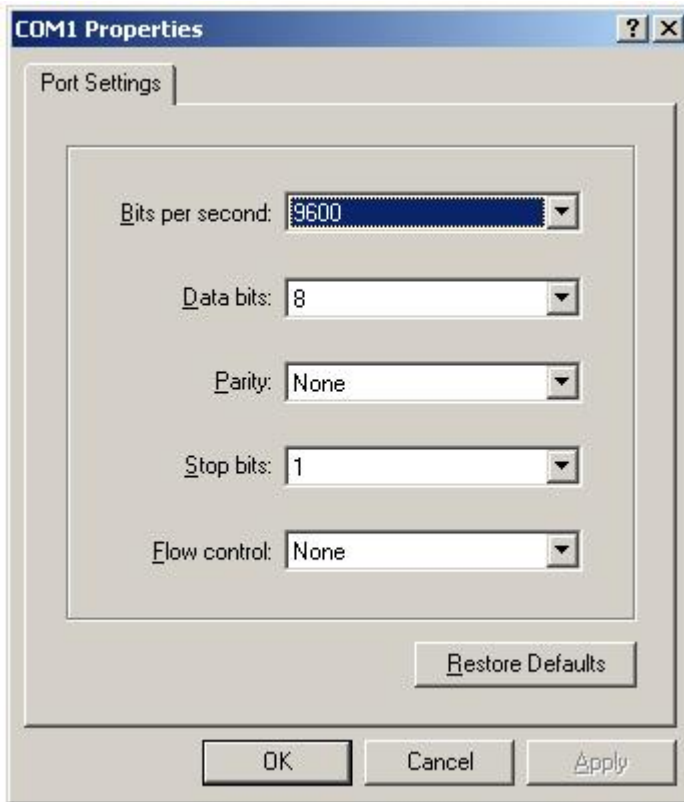
## **Syringe Pump Controller Program** (Freeware not supported)

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